# **Application Note AFS-101**

## **Determination of Mercury in Water Samples by HG-AFS**

Mercury and its compounds are highly toxic materials and can be transferred into the human body via water; the concentration of mercury in water generally is very low. With the superior sensitivity of atomic fluorescence spectrometer, direct measurement of mercury in water sample is possible.

## 1. Major equipment and reagents.

## Mercury stabilization solutions:

Dissolve 0.5g K<sub>2</sub>Cr<sub>2</sub>O<sub>7</sub> in 950 mL water, add 50ml nitric acid.

#### Mercury standard solutions:

Prepare or commercial available mercury standard at 100ppm, diluted to 10ppb with mercury stabilization solution as above.

KMnO₄ solution, 50g/L.

Potassium persulfate solution, 50 g/L

#### SnCl, solution:

Add 100g SnCl2 in 100 mL concentrated HCl, water bath to totally dissolved, add water to 1000 mL.

NaCl-NH<sub>2</sub>OH solution, 120 g/L

AI3300 atomic fluorescence spectrometer with Hg lamp.

High pure distilled water.

High pure argon (>99.99%)

#### 2. Method

Place 25mL water sample in a 50mL flask, add 4mL concentrated  $H_2SO_4$ , 2.5mL KMnO $_4$  solution, and 1.0mL potassium persulfate solution, in a boiling water bath for 1 hr. Take the flask from the bath and cool it down. Before measurement, add NaCl-NH $_2$ OH solution to the color of KMnO $_4$  until the color disappears, then add water until the total volume is 50mL. Use the final solution for measurement.

### 3. Instrument parameters

Carrier gas 300mL/min
Shield gas 800mL/min
HCL current 30mA
PMT voltage 360V
Integration time 6 s
Pump speed 40 r/min

Reducing reagent solution 10% SnCl<sub>2</sub> in 10% HCl (V/V)

#### 4. Results

This method gives: Detection limit: 0.25ppb, Recovery rate: 98~117%

Relative standard deviation: 2~6%

